(90 × 210mm Size)

Code No. 293-77601 (10 purifications)
299-77603 (2 purifications)

For Genetic Research
MagCapture™ Exosome Isolation Kit PS

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Exosomes and other extracellular vesicles (EVs) are small membrane vesicles containing protein, mRNA, microRNA, DNA, and lipids, which are secreted by various cells and are stable in body fluids including blood, saliva, urine, cerebrospinal fluid (CSF) and breast milk. These EVs have been recognized as messengers of cell-to-cell communication, and biomarkers for various diseases. Conventionally, ultracentrifugation, affinity purification using antibody to surface antigen, and precipitation with polymer reagents are used for isolation of exosomes. However, these methods are not fully satisfying in all performances criteria such as recovery efficiency, purity and operability.

MagCapture™ Exosome Isolation Kit PS adopts a novel affinity purification method using magnetic beads and phosphatidylserine (PS)-binding protein (PS affinity method). This kit can easily isolate high purity exosomes and other EVs from cell culture medium and body fluids at high yield by a normal microcentrifuge. If higher purity exosomes are needed, please use the 10,000 × g supernatant as sample. This kit enables the isolation of exosomes and other EVs as intact forms because the captured EVs are eluted from magnetic beads with metal-chelating reagent at neutral pH. The isolated intact exosomes and other EVs can be used for various applications including electron microscopic analysis, nanoparticle tracking analysis, administration of EVs and analysis of molecular constituents such as proteins, lipids, or nucleic acids.

【Features】

- High purity exosomes and other EVs are isolated by a novel affinity purification method (PS affinity method).
- The purity and yield of exosomes are higher than ultracentrifugation.
- This kit enables the purification of exosomes and other EVs as intact forms, and can be used for various applications.
- Exosomes can be isolated with magnetic beads.
- No ultracentrifugation is required.

【Kit contents】

This kit includes 7 components.

· 2 purifications

(1) Streptavidin Magnetic Beads	120 μL × 1 tube
(2) Biotin-labeled Exosome Capture	20 μL × 1 tube
(3) Exosome Capture Immobilizing Buffer	7 mL × 1 bottle
(4) Exosome Binding Enhancer (× 500)	100 μL × 1 tube
(5) Washing Buffer	30 mL × 1 bottle
(6) Exosome Elution Buffer	1 mL × 1 bottle
(7) Reaction Tubes	4 tubes

– 1/20 –

· 10 purifications

(1) Streptavidin Magnetic Beads	600 μL × 1 tube
(2) Biotin-labeled Exosome Capture	100 μL × 1 tube
(3) Exosome Capture Immobilizing Buffer	35 mL × 1 bottle
(4) Exosome Binding Enhancer (× 500)	500 μL × 1 tube
(5) Washing Buffer	75 mL × 2 bottles
(6) Exosome Elution Buffer	5 mL × 1 bottle
(7) Reaction Tubes	22 tubes

【Storage】

Store at 2-10 °C

【Additional required materials】

Reagents :

- 1) TBS (As necessary)

Equipment :

- 1) Microcentrifuge (Max > 10,000 × g)
- 2) Vortex mixer
- 3) Tabletop centrifuge
- 4) Magnetic stand
- 5) Rotator or tube mixer
- 6) Micro pipette
- 7) Pipette tip
- 8) Microcentrifuge tube (1.5 mL)
- 9) Centrifuge tube (15 mL) (As necessary)
- 10) Centrifuge tube (50 mL) (As necessary)
- 11) Ultrafiltration unit (Vivaspin6, M. W. cut off : 100K, Code No. VS0641 or Vivaspin20, M. W. cut off : 100K, Code No. VS2041) (As necessary)

【Precaution for use】

1. Equipment

Please use sterile or DNase and RNase free microcentrifuge tubes and pipette tips. We recommend the use of gloves and a mask to avoid contamination of DNase and RNase.

2. Reagents


If you analyze the nucleic acid (RNA, DNA) after isolation of exosomes and other EVs, please handle reagents carefully and keep them as sterile as possible. Additionally, if you analyze the mass spectrometry of proteins after isolation of exosomes and other EVs, please be careful of contamination from other proteins.

3. Handling of biohazardous waste

Please treat human serum, human plasma, and human tissue samples as infectious samples. Biohazardous wastes, which include waste liquid and equipment such as microcentrifuge tubes, pipette tips, and gloves, must be disposed of according to the guidelines of the institution.

– 2/20 –

【Outline of procedure】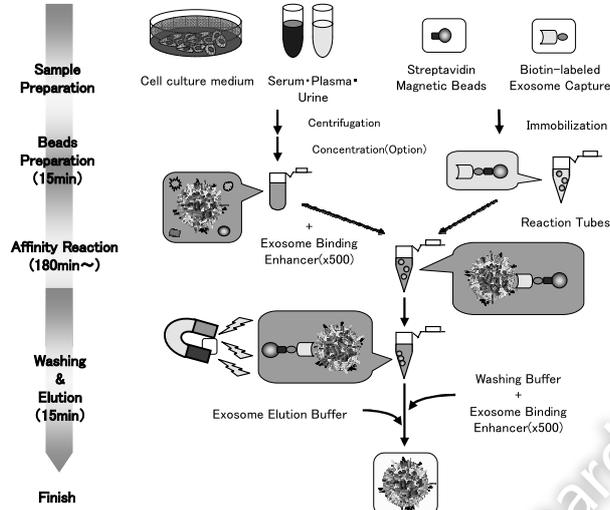

【Procedure】

1. Preparation of sample

This is the section to prepare samples. When exosomes and other large EVs are needed, prepare **1,200 × g supernatant** as a sample. Additionally, when highly purified exosomes are needed, prepare **10,000 × g supernatant** as a sample. This protocol for sample preparation is set for cell culture medium, serum, and plasma. When other body fluids are used, please examine the appropriate pretreatment protocol by referring to the protocol for serum and plasma.

In the case of cell culture medium

- 1) Culture the cells under the appropriate condition^{※1}.
- 2) Harvest cell culture medium.
- 3) Centrifuge the cell culture medium for 5 minutes at 300 × g and 4 °C to remove cells and debris.
- 4) Transfer the supernatant from step 3) into a new tube.
- 5) Centrifuge 4) at 1,200 × g for 20 minutes at 4 °C to remove the cell debris.
- 6) Transfer the supernatant from step 5) into a new tube. (**1,200 × g supernatant**)
- 7) Centrifuge 6) at 10,000 × g for 30 minutes at 4 °C to remove the large EVs.
- 8) Transfer the supernatant from step 7) into a new tube. (**10,000 × g supernatant**)

【Optional】

If the volume of **1,200 × g supernatant** or **10,000 × g supernatant** is over 1 mL, concentrating the sample by using ultrafiltration unit (Vivaspin6, M. W. cut off: 100K, Code No. VS0641 or Vivaspin20, M. W. cut off: 100K, Code No. VS2041) is recommended. Please concentrate until the sample is less than or equal to 1 mL in accordance with the manufacturer's instruction manual. The maximum sample volume of ultrafiltration treatment is 50 mL (50-fold concentration). Exosomes and other EVs are efficiently recovered from concentrated sample.

※1 Please refer to appropriate culture conditions depending on cell lines. To ensure that isolated exosomes which originate from your cells, please use exosome-depleted FBS. Exosome-depleted FBS is prepared by ultracentrifugation (Example: at 110,000 × g for 18 hours) or is commercially available.

In the case of serum

- 1) Centrifuge the serum at 1,200 × g for 20 minutes at 4 °C to remove the cell debris.
- 2) Transfer the supernatant from step 1) into a new tube. (**1,200 × g supernatant**^{※2})
- 3) Centrifuge 2) at 10,000 × g for 30 minutes at 4 °C to remove the large EVs.
- 4) Transfer the supernatant from step 3) into a new tube. (**10,000 × g supernatant**^{※2})

※2 In the case of serum, 1,200 × g supernatant and 10,000 × g supernatant are unsuitable for concentration with ultrafiltration unit.

In the case of plasma

Please prepare as much Washing Buffer as plasma.


- 1) Centrifuge the plasma at 1,200 × g for 20 minutes at 4 °C to remove the cell debris.
- 2) Transfer the supernatant from step 1) into a new tube. (**1,200 × g supernatant**^{※3})
- 3) Centrifuge 2) at 10,000 × g for 30 minutes at 4 °C to remove the large EVs.
- 4) Transfer the supernatant from step 3) into a new tube. (**10,000 × g supernatant**^{※3})
- 5) Add 1:100 volume of Exosome Binding Enhancer (× 500) to Washing Buffer^{※4}.
- 6) Add the same amount of Washing Buffer including Exosome Binding Enhancer (× 500) into a plasma sample from step 2) or step 4).
- 7) Incubate with rotator for 10 minutes at room temperature. ... **A sample including Exosome Binding Enhancer (× 500)**
- 8) Proceed to step 2) of section 3.

※3 In the case of plasma, 1,200 × g supernatant and 10,000 × g supernatant are unsuitable for concentration with ultrafiltration unit.

※4 Please adjust the volume of Washing Buffer including Exosome Binding Enhancer (× 500) depending on the volume of plasma sample.

【Flowchart of Sample Preparation (Section 1)】

Cell Culture Medium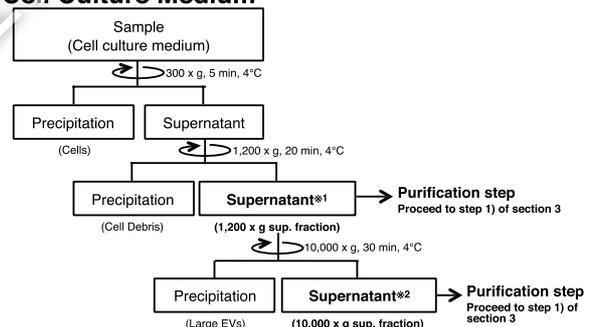

Concentration (Optional)^{※3}


~50 mL (1,200 × g Sup. fraction or 10,000 × g Sup. fraction)

Ultrafiltration and concentration (VivaSpin20, M. W. cut off: 100K)


1 mL → Purification step

(~50-fold concentrated)

Serum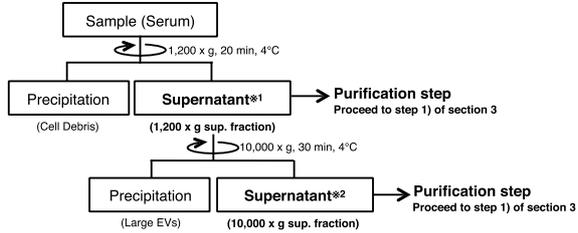

Plasma

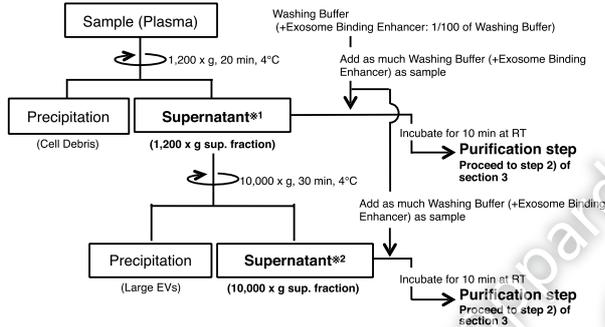※1 When **exosomes and large EVs** are needed, use 1,200 × g sup. fraction as sample.

※2 When **exosomes** are needed, use 10,000 × g sup. fraction as sample.

※3 This is optional protocol for large-scale volume (~50 mL) of cell culture medium.

2. Immobilization of Exosome Capture


This is the section to immobilize Biotin-labeled Exosome Capture reagent to Streptavidin Magnetic Beads. Please use the 1.5 mL Reaction Tube (included) in the downstream experiment which requires magnetic beads operation※5.

- 1) Transfer 60 μL ※6 of Streptavidin Magnetic Beads into a new 1.5 mL Reaction Tube.
- 2) Place the 1.5 mL Reaction Tube on the magnetic stand for 1 minute to collect the beads. Remove the supernatant after separation of magnetic beads and supernatant.
- 3) Add 500 μL of Exosome Capture Immobilizing Buffer into the 1.5 mL Reaction Tube from step 2) and suspend it by vortexing.
- 4) Spin down 3), then place it on the magnetic stand for 1 minute. Remove the supernatant after separation of magnetic beads and supernatant.
- 5) Add 500 μL of Exosome Capture Immobilizing Buffer and 10 μL of Biotin-labeled Exosome Capture into the 1.5 mL Reaction Tube from step 4) and mix by vortexing.
- 6) Mix for 10 minutes at 2-10 $^{\circ}\text{C}$ with rotator or tube mixer.
- 7) Spin down 6), then place it on the magnetic stand for 1 minute. Remove the supernatant after separation of magnetic beads and supernatant.

- 8) Add 500 μL of Exosome Capture Immobilizing Buffer into the 1.5 mL Reaction Tube from step 7) and mix by vortexing.
- 9) Spin down 8), then place it on the magnetic stand for 1 minute. Remove the supernatant after separation of magnetic beads and supernatant.
- 10) Repeat steps 8)-9) twice. . . . **Exosome Capture-immobilized beads**

- ※5 Please only use the 1.5 mL Reaction Tube included in this kit because only these tubes can generate the high recovery rate of extracellular vesicles by preventing non-specific adsorption of magnetic beads to the tube wall.
- ※6 Please adequately suspend the magnetic beads by vortexing. Make sure homogeneous suspension of the magnetic beads is achieved before use.

[Flowchart of immobilization of Exosome Capture (Section 2)]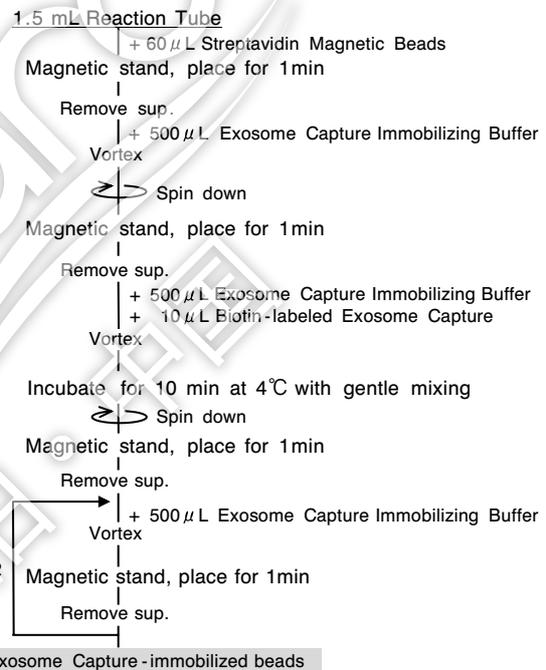

3. Affinity Reaction

This is the affinity reaction section of the Exosome Capture-immobilized beads with exosomes and other EVs in sample. This protocol is based on the use of 1.5 mL Reaction Tube (max volume : 1 mL). When exosomes and large EVs are needed, use **1,200 × g supernatant** as sample. Additionally, when highly purified exosomes are needed, use **10,000 × g supernatant** as sample.

- 1) Transfer sample (max volume : 1 mL)※7 to a new 1.5 mL microcentrifuge tube (not included in this kit) and add 1:500 volume of Exosome Binding Enhancer (× 500) to sample. Mix by vortexing. . . . **A sample including Exosome Binding Enhancer (× 500)**

- 2) Spin down **A sample including Exosome Binding Enhancer (×500)**, then transfer it into the 1.5 mL Reaction Tube containing **Exosome Capture-immobilized beads**. Mix by vortexing.
- 3) Mix for 3 hours~ at 2-10 °C with rotator or tube mixer.
- 4) Spin down 3), then places it on the magnetic stand for 1 minute. Remove the supernatant^{※8} after separation of magnetic beads and supernatant. . . . **EVs-binding beads**

[Optional]

When the sample volume is over 1 mL, please refer to the following protocol to use 15 mL centrifuge tube (max volume: 10 mL). However, the increase of sample volume may decrease recovery rate. Therefore, we recommend the use of the Exosome Capture-immobilized beads repeatedly in order to recover extracellular vesicles from used sample (refer to ※16). Please repeat sections 3-5 (up to 4 times). Additionally, in the case of cell culture medium, we recommend a concentration less than or equal to 1 mL by using ultrafiltration unit in accordance with above-described protocol.

- 1) Transfer sample (max volume: 10 mL) to a new 15 mL centrifuge tube^{※9} (not included in this kit) and add 1:500 volume of Exosome Binding Enhancer (×500) to sample. Mix by vortexing. . . . **A sample including Exosome Binding Enhancer (×500)**
- 2) Transfer 1 mL of **A sample including Exosome Binding Enhancer (×500)** into the 1.5 mL Reaction Tube^{※10} containing **Exosome Capture-immobilized beads**. After suspending the beads, transfer all of the beads in suspension back to the 15 mL centrifuge tube. Mix by vortexing.
- 3) Mix for 3 hours~ at 2-10 °C with rotator or tube mixer.
- 4) Centrifuge 3) at 1,000 × g for 3 minutes at 4 °C.
- 5) Remove the supernatant^{※8} carefully and leave 1ml of the supernatant and bead pellets.
- 6) After suspending sample from step 5), transfer all of the beads in suspension into the 1.5 mL Reaction Tube from 2).
- 7) Spin down 6), then place it on the magnetic stand for 1 minute to collect the beads. Remove the supernatant^{※8} after separation of magnetic beads and supernatant. . . . **EVs-binding beads**

※7 Add the appropriate volume of TBS buffer to each sample to reach a volume of 0.5 mL to obtain the better mixture of the Exosome Capture-immobilized beads with sample. (Example : 100-200 μL→500 μL)

※8 When further recovery of remaining extracellular vesicles is needed, please store the used supernatant.

※9 Please use a 15 mL centrifuge tube in which the non-specific adsorption of magnetic beads to the tube wall is low. (Example : Code No. 430791, Corning)

※10 Please store this 1.5 mL Reaction Tube after transferring the beads into the 15 mL centrifuge tube, because it is used again at the optional step 6) of section 3.

4. Washing of EVs-binding beads

This is the washing section of **EVs-binding beads**. Please use Washing Buffer after adding Exosome Binding Enhancer (×500) certainly^{※11}.

- 1) Add 1:500 volume of Exosome Binding Enhancer (×500) to the required Washing Buffer^{※12} and mix by vortexing^{※13}.
- 2) Add 1 mL of Washing Buffer (+ Exosome Binding Enhancer (×500)) into the 1.5 mL Reaction Tube containing **EVs-binding beads** and mix by vortexing.
- 3) Spin down 2), then places it on the magnetic stand for 1 minute. Remove the supernatant after separation of magnetic beads and supernatant.
- 4) Repeat steps 2)-3) twice.
- 5) Spin down again the 1.5 mL Reaction Tube from step 4), then places it on the magnetic stand for 1 minute. Remove the supernatant completely after separation of magnetic beads and supernatant^{※14}. . . . **Washed EVs-binding beads**

※11 If the addition of Exosome Binding Enhancer (×500) to Washing buffer is omitted, captured exosome and EVs will be eluted from the magnetic beads by Washing Buffer during washing section.

※12 The required volume of Washing Buffer is 3 mL per reaction.

※13 Please prepare Washing Buffer (+ Exosome Binding Enhancer (×500)) just before use, because constituents may be precipitated after the addition of Exosome Binding Enhancer (×500) overtime.

※14 Please remove Washing Buffer (+ Exosome Binding Enhancer (×500)) in the tube completely, because remaining buffer causes reduction of elution efficiency of captured exosomes and other EVs.

5. Elution of extracellular vesicles

This is the elution section of extracellular vesicles from **Washed EVs-binding beads**.


- 1) Add 50 μL of Exosome Elution Buffer into the 1.5 mL Reaction Tube containing **Washed EVs-binding beads** and mix by vortexing^{※15}.
- 2) Spin down 1), then places it on the magnetic stand for 1 minute. Transfer the supernatant into a new 1.5 mL microcentrifuge tube (not included in this kit) after separation of magnetic beads and supernatant.
- 3) Add 50 μL of Exosome Elution Buffer into the 1.5 mL Reaction Tube from step 2) again, and mix by vortexing^{※15}.
- 4) Spin down 3), then places it on the magnetic stand for 1 minute. Transfer the supernatant into the 1.5 mL microcentrifuge tube from step 2) after separation of magnetic beads and supernatant, and then pool together^{※16, 17}. (Total: 100 μL)

[Optional]

In the case of RNA purification from eluted extracellular vesicles, please add 100 μL of Exosome Elution Buffer to pooled eluate to prepare 200 μL of sample in total. Then, extract RNAs from extracellular vesicles with Wako's microRNA Extractor SP Kit (Code No. 295-71701) in accordance with the instruction manual.

- ※15 When the aggregates of magnetic beads are remaining, please suspend magnetic beads by tapping and/or pipetting until it reaches a uniform consistency.
- ※16 **Exosome Capture-immobilized beads** are reusable after eluting EVs (up to 4 times). Therefore, when you need to recover more remaining EVs from samples, please repeat the section 3-5 to increase the recovered amount. After recovering the eluate, please add the stored supernatant of section 3 (refer to ※8) into the 1.5 mL Reaction Tube containing **Exosome Capture-immobilized beads** from step 4) of section 5 and repeat the section 3-5 as necessary. This kit includes the required reagents to reuse a maximum of 4 times.
- ※17 There is a possibility that the pooled eluate have a small amount of magnetic beads. When extracellular vesicles are analyzed by Nanoparticle Tracking Analysis (NTA) or electron microscopic analysis, please filter pooled eluate by using 0.45 μm pore size filter unit (Example: Ultrafree[®] MC 0.45 μm, Code No. UFC30HV25, Millipore[®]) before analysis as needed.

[Flowchart of Affinity Purification (Section 3-5)]

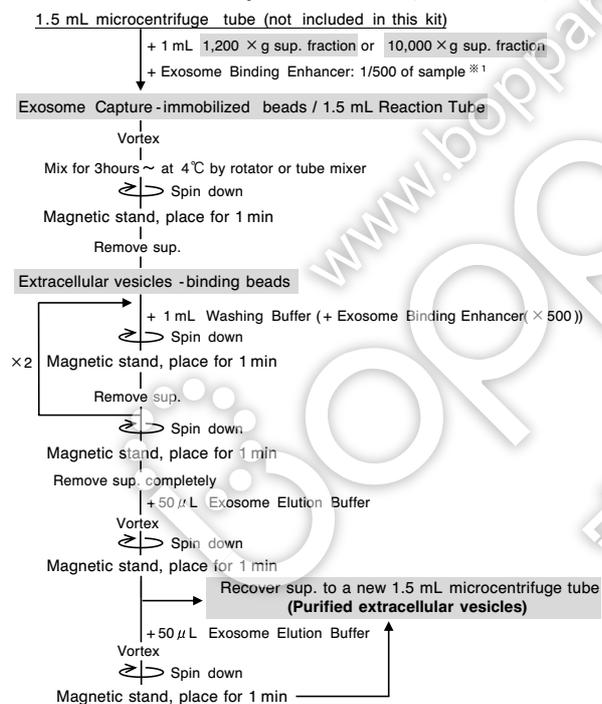※1 In the case of plasma, no need to add the Enhancer again.

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Code No.	Description	Size
295-71701	microRNA Extractor SP Kit	50 reactions
317-90175	10 × TBS (pH 7.4)	500 mL

【Reference】

W. Nakai, *et al.*, *Sci Rep* **6**, 33935 (2016).

Wako Pure Chemical Industries, Ltd.

1-2, Doshomachi 3-Chome, Chuo-Ku, Osaka 540-8605, Japan
 Telephone : + 81-6-6203-3741
 Facsimile : + 81-6-6201-5964
<http://www.wako-chem.co.jp>

Wako Chemicals USA, Inc.

1600 Bellwood Road
 Richmond, VA 23237
 U.S.A.
 Telephone : + 1-804-271-7677
 Facsimile : + 1-804-271-7791
<http://www.wakousa.com>

Wako Chemicals GmbH

Fuggerstrasse 12
 D-41468 Neuss
 Germany
 Telephone : + 49-2131-311-0
 Facsimile : + 49-2131-311100
<http://www.wako-chemicals.de>

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MagCapture™ Exosome Isolation Kit PS

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