

IHC Protocol (Paraffin Section) Anti Iba1, Rabbit

100% xylene, 5 min×2
100% Ethanol, 5 min ×2
0.3% H₂O₂ in Methanol, 20 min, (inactivation of endogenous peroxidase)
90% Ethanol , 2 min
70% Ethanol , 2 min
H₂O 3 min
PBS 5 min×2
↓
Citrate Buffer (pH6.0) 90°C, 9 min
Incubate to RT
↓
PBS 5 min×2
↓
2% BSA in PBS RT, 60min (blocking)
↓
PBS 5 min×2
↓
Anti Iba1, Rabbit (1:1,000) O/N
↓
PBS 5 min×2
↓
Secondary antibody (Biotinylated anti Rabbit IgG), 60 min, RT
↓
PBS 5 min×2
↓
ABC, 1 hour, RT
↓
PBS, 5 min×2
↓
DAB solution
↓
PBS, 5 min×2
↓
Mount

产品列表

产品编号	产品名称	包装
019-19741	小胶质细胞/巨噬细胞特异性蛋白抗体（免疫组化） Anti Iba1, Rabbit (for Immunocytochemistry)	50 µg

